



Guide to trapping moulds / fungi - using Petri dishes (agar plates)

- Obtain the required number of Petri dishes containing nutrient agar from PLANTwise
- Place plates throughout the building at sites where moulds are suspected to be air-borne - LABEL with location and date
- Last thing at night, remove the Petri dish lids
- First thing in the morning, replace the lids and secure the lids with adhesive tape (2 -3 short strips from lid to base)
- Return the plates to the PLANTwise (by courier or delivery) for incubation, enumeration and mould identification

Details:

Your reference:

Date:

Location:

.....

.....

Contact details:

Collector:

Contact details:

.....

.....

.....

Email address:

.....

Notes:

.....

.....

.....

For Lab use:

Our Reference:

Date received:

SEND TO:

PLANTwise Services Ltd

PO Box 69181

Lincoln 7640

NEW ZEALAND

Work: 03 3257 240

Mobil: 027 2227 435

COURIER TO:

55a Lochee Rd, Christchurch 8041

Email: harveyi@plantwise.co.nzWeb: www.plantwise.co.nz